

Ortho-Jet™ BCA

Product Identification

Self Curing Acrylic Resin

Package Contents

REF B1320	Net Wt. 100g
REF B1330	Net Wt. 454g
REF B1350	Net Wt. 2.3kg
REF B1303	Net Vol. 118ml
REF B1304	Net Vol. 236ml
REF B1306	Net Vol. 946ml
REF B1323	Net Wt. 100g; Net Vol. 118ml; glass dropper
REF B1334	Net Wt. 454g; Net Vol. 236ml; glass dropper; dispenser bottle
REF B1356	Net Wt. 2.3kg; Net Vol. 946ml

Intended Purpose

This product is intended for research use in potting, embedding and cementation.

Storage Conditions/Instructions

Close container tightly after use. Store in original container, in a cool, dry, well-ventilated location, away from sources of ignition and light.



F - Highly Flammable

Operating/Handling Instructions

Liquid may cause skin irritation. Avoid contact with skin, eyes and clothing. Wear protective gloves, safety glasses with side shields or chemical splash goggles. Avoid prolonged breathing of vapors. Use with adequate ventilation. Do not smoke in the presence of open containers. See Material Safety Data Sheet for complete information. MSDS is available at www.langdental.com or by contacting manufacturer.



Xn - Harmful

Warning: flammable liquid. Contains: Methyl Methacrylate Monomer, stabilized. Flash Point: 52.7°F (11.5°C), NFPA, NPCA-HMIS #: H-2,F-3,R-2.

Other Materials Needed

1. Mixing cup
2. Stirring instrument

Instructions for Use

For consistent results always shake liquid before use.

1. Mix 2 parts by volume of Ortho-Jet BCA powder with 1 part by volume of Ortho-Jet BCA liquid. (For lower viscosity, mix 3 parts by volume of powder with 2 parts liquid).
2. Blend mixture thoroughly for 15-20 seconds.
3. Apply mixture as needed.
4. Allow to cure at room temperature.

Disposal

Dispose of unused material in accordance with local, state, and national legislation.

Warranty

Manufacturer warrants this product to be free from manufacturing defects for one year from date of purchase. Warranty is void if product is not used for intended purposes, not used as instructed or not stored in original container.

Lang Dental Mfg. Co., Inc.
Wheeling, IL 60090-0969
www.LangDental.com
Made in USA
Ph 847-215-6622
Fax 847-215-6678

The logo for Lang Dental features the word "LANG" in a bold, italicized, sans-serif font. The letters are contained within a horizontal rectangular frame. A diagonal line extends from the bottom-left corner of the frame, passing through the letter 'A' and continuing downwards and to the left.