

Ortho-Jet™ BCA

Intended Purpose

Potting, embedding and cementation for research

Other Materials / Equipment Needed

Mixing cup
Stirring instrument

Instructions for Use



For consistent results, always shake liquid before use, as catalyst may have separated from monomer.

1. Mix 2 parts by volume of Ortho-Jet™ BCA powder with 1 part by volume of Ortho-Jet™ BCA liquid. (For lower viscosity, mix 3 parts by volume of powder with 2 parts liquid).
2. Blend mixture thoroughly for 15-20 seconds.
3. Apply mixture as needed.
4. Allow to cure at room temperature.

Handling Precautions

Contains methyl methacrylate

Danger. Highly flammable liquid and vapor. Keep away from heat/sparks/open flames/hot surface. - No smoking.

Causes skin irritation. May cause an allergic skin reaction. May cause respiratory irritation.

Wear protective gloves/protective clothing/eye protection/ face protection.

Keep material in original container tightly closed in upright position. Store in a well-ventilated place. Keep cool.

Disposal

Dispose of contents and container in accordance with local, regional and national regulations.


Warranty

Manufacturer warrants this product to be free from manufacturing defects for one year from date of purchase only from an authorized reseller. Warranty is void if product is not used for intended purpose, not used as instructed or not stored in original container.



Lang Dental Mfg. Co., Inc.
175 Messner Dr
Wheeling, IL 60090
USA

Ph +1 847-215-6622
Fax +1 847-215-6678
www.LangDental.com

Made in USA 

© 2023 Lang Dental Mfg. Co., Inc. LANG is the registered trademark of Lang Dental Mfg. Co., Inc. ORTHO-JET is the trademark of Lang Dental Mfg. Co., Inc. All rights reserved.

LDD B13 R. 1 2023-01